



SETUP AND USE OF VP 710J-AIR (AIR MOTOR TUMBLE STIRRER)

Setup:

1. Connect a hose from port labeled "A" on the tumble stirrer to the control unit lubricator port labeled "A".
2. Connect a hose from port labeled "B" on the tumble stirrer to the control unit speed control port labeled "B".
3. Connect the hose with the industrial-shape hose coupling to the control unit port labeled "AIR SOURCE", the coupling on the other end of the hose is to be connected to a compressed air source.
4. Fill about 75% of the lubricator with pneumatic oil.

Use:

1. Turn the main valve ON in the control unit
2. Adjust the pressure regulator the desired pressure use (<150 psi).
3. Adjust the speed using the speed control (notice label at the bottom of speed control valve)
4. To improve speed, try to keep the air muffler as close as possible to the exhaust port.