



**IN THE RARE EVENT THAT AN E-CLIP FALLS OFF,  
HERE ARE SOME STEPS TO ATTACH IT BACK ON.**

1. Grab the E-clip with needle nose pliers.
  - a. The jaws of the pliers should be just at the edge of the E-clip, leaving the center open.
  - b. Position the E-clip such, so that the shiny somewhat rounded side of the E-clip is facing up.
2. Line up the groove on Pin with the slot of the E-clip.
3. Press the Pin into the E-clip so the groove slides inside the opening of the E-clip.
4. Make the sure E-clip is secure and straight in the groove.